# MVE Operating Instruction Manual for Liquid Nitrogen Dewars

<b>Table of Contents</b>	<b>Page</b>
English	2-5
English (SI Version)	6-9
German	10-13
Spanish	14-17
Italian	18-21
French	22-25

**Chart Bio-Medical Division,  
Chart Industries, Inc.**  
1800 Sandy Plains Industrial Parkway  
Suite 316  
Marietta, GA 30066-6363 USA  
[www.chartbiomed.com](http://www.chartbiomed.com)

**Chart Bio-Medical Products,**  
112 Windmill Road  
Sunbury on Thames  
Middlesex TW167HB  
United Kingdom

**Chart Australia Pty Ltd.  
Sidney Business & Technology Centre**  
Unit 43 / 2 Railway Parade  
Lidcombe NSW 2141, Australia

**M.D.D. Representative:**  
Medical Product Services  
Borngasse 20  
35619 Braunfels, Germany

Ref 11624417 Rev D 6/05

**GENERAL DESCRIPTION**

The SC, XC and LAB series container is a vacuum insulated container of aluminum with fiberglass neck construction providing the highest efficiency possible in liquid nitrogen storage. Use the container for inert fluids only. Liquid oxygen is not compatible with fiberglass material and must not be stored. A sharp blow to the outer vessel can damage the neck tube or start a vacuum leak. Use caution and common sense in handling the container. The life expectancy of the CryoSystem Series is five (5) years. SC, XC, and Lab Series have a life expectancy of ten (10) years.

Upon receipt of the container, examine both the container and packaging for any evidence of damage during shipping. Contact the shipping agent if any damage is apparent. Watch after the first fill for any signs of vacuum loss, such as frost or sweating on the outside jacket. (Some frost near the tip just after filling is normal).

**A. SAFETY**

-  To avoid injury by frostbite, use extreme care whenever handling liquid nitrogen, liquid nitrogen storage, transfer vessels or any objects which have come in contact with liquid nitrogen.
  - Leave no areas of skin exposed.
  - Always wear proper safety attire over clothing: face shield, cryogenic gloves, cryogenic apron.
  - Never overfill liquid nitrogen vessels.
-  Always keep liquid nitrogen vessel in an upright position.
-  Do not tightly seal liquid nitrogen container or prevent nitrogen gas from escaping.
-  Use extreme care to prevent spilling and splashing liquid nitrogen during transfer.
- Immediately remove any clothing or safety attire on which liquid nitrogen has been spilled.
- Get immediate medical attention for any frostbite injuries due to liquid nitrogen.

 **WARNING:** The venting of nitrogen vapors will deplete oxygen in the air, possibly leading to asphyxiation or even death. **DO NOT** store or use liquid container in areas that have poor ventilation.

**B. FILLING INSTRUCTIONS**

- To avoid damage to the aluminum cryogenic vessel which may result in premature vacuum loss, it is important to follow this procedure during the addition of liquid nitrogen to a warm vessel and on subsequent additions.
1. Slowly pour liquid nitrogen into new or warm vessels with a funnel or transfer line when possible.
  2. Allow liquid nitrogen to settle in vessel for 2 hours to completely cool inner.
  3. Fill the vessel to the desired level after the 2 hr settling (cooldown) time.
  4. If filling the dewar from a pressurized source, make sure that the source tank is at a low pressure (22 psi or below).
  5. If transfer hose is used for extracting liquid nitrogen from a pressurized liquid source always use a phase separator on the end of the hose.
  6. Remember to always wear proper safety attire over clothing; face shield, cryogenic gloves, and apron.
  7. Never overfill your dewar with liquid nitrogen. Overfilling the tank may cause pumpout components to leak leading to immediate or premature vacuum failure to occur.

**C. MEASURING LIQUID NITROGEN QUANTITY**

1. Use wooden or plastic dipstick. Never use a hollow tube to measure liquid nitrogen.
2. Level will be indicated by frostline, which develops when dipstick is removed.

**D. LIQUID WITHDRAWAL**

1. Liquid withdrawal from LAB units must be accomplished by pouring or utilizing a withdrawal device. Withdrawal device pressurizes to approximately 5 psi and the pressure forces liquid up the withdrawal tube out the valve.
2. Always wear proper safety attire; shield, gloves and apron.

**WARRANTY ACTIVATION**

**IMPORTANT! PLEASE FILL OUT THIS REGISTRATION CARD TO ACTIVATE YOUR WARRANTY**

YOUR NAME \_\_\_\_\_  
 ADDRESS \_\_\_\_\_  
 CITY \_\_\_\_\_ STATE \_\_\_\_\_ ZIP \_\_\_\_\_  
 NAME OF DISTRIBUTOR \_\_\_\_\_  
 ADDRESS \_\_\_\_\_  
 SERIAL NUMBER \_\_\_\_\_ DATE PURCHASED \_\_\_\_\_

*What is being stored in the container?*

- Semen                       Embryos                       Skin                       Vaccines                       Blood  
 Bone Marrow                       Cell Lines                       Other

*Was the new container purchased for*

- Replacement                       New Lab                       Project                       Additional Storage

*How did you become aware of MVE products?*

- Advertisement                       Sales Distributor                       Colleague                       Trade Show  
 Thomas Register                       Internet                       Other \_\_\_\_\_

*What Journal/Magazines are you currently receiving?*

(Rate 1 as most useful, 3 as less useful)

1 \_\_\_\_\_ 2 \_\_\_\_\_ 3 \_\_\_\_\_

*How important were the following features that lead to the purchase of this product?*

	Very Important							Not Important		
Quality of freezer	10	9	8	7	6	5	4	3	2	1
MVE reputation	10	9	8	7	6	5	4	3	2	1
Price	10	9	8	7	6	5	4	3	2	1
Storage Capacity	10	9	8	7	6	5	4	3	2	1
Warranty	10	9	8	7	6	5	4	3	2	1
Standard Features	10	9	8	7	6	5	4	3	2	1
Customer Service	10	9	8	7	6	5	4	3	2	1
Distributor Reputation	10	9	8	7	6	5	4	3	2	1
Colleague Referral	10	9	8	7	6	5	4	3	2	1
Temperature Under Lid	10	9	8	7	6	5	4	3	2	1
Nitrogen Consumption	10	9	8	7	6	5	4	3	2	1
Serviceability	10	9	8	7	6	5	4	3	2	1
Other	10	9	8	7	6	5	4	3	2	1

*How satisfied are you with your local MVE representative?*

	Very Satisfied							Not Satisfied		
	10	9	8	7	6	5	4	3	2	1

Comments \_\_\_\_\_

*Overall, how satisfied are you with your MVE container?*

	10	9	8	7	6	5	4	3	2	1

Comments \_\_\_\_\_

Please make any additional comments or suggestions \_\_\_\_\_

**CHART BIO-MEDICAL PRODUCTS**  
**ATTN: CUSTOMER SERVICE MANAGER**  
**1800 Sandy Plains Industrial Parkway**  
**Suite 316**  
**Marietta, GA 30066-6363**

PLACE  
STAMP  
HERE

---

---

---

---

*Thank You!* You have just purchased the finest cryobiological storage container manufactured today. With Chart Bio-Medical Division's experience and reputation, you can be assured of the most advanced technology and unsurpassed quality throughout our product. These containers have been designed, manufactured and tested to achieve maximum efficiency, along with durability to give many years of trouble-free service. If you have any questions regarding this product please contact Chart Bio-Medical.

## WARRANTY

MVE warrants to the original purchaser ("Purchaser") that each new Liquid Nitrogen Dewar (each a "Product"), from the date of shipment, shall be free from defects in materials and workmanship for a period of two (2) years, except as provided below, and dewar vacuum integrity shall be free from defects in materials and workmanship as listed below, except as provided below:

Spectrum and CryoSystem Series, ET11, ET24, A-200, Workhorse 2003: 3 years / XC, SC, and Lab Series, ET 23-JLM: 5 years

MVE warrants all repaired equipment and factory blemished equipment shall be free of defects in materials and workmanship for a period of (90) days from the date of shipment and warrants all dewar vacuum integrity shall be free from defects in materials and workmanship for a period of (1) year from date of shipment. All accessory items are warranted for (90) days from date of shipment to be free of defects in material and workmanship. Purchaser agrees that before this limited warranty shall become effective, Purchaser shall fully inspect each Product within ten (10) days of delivery and before such Product is put to use. Purchaser also agrees to operate the Product in accordance with MVE's operating instructions and that failure to do so shall void this limited warranty. Purchaser further agrees that any claim for breach of warranty must be made in writing within 60 days of discovery of a purported defect. MVE will not be responsible for any alleged breach of warranty, which, as a result of MVE's inspection, MVE determines to have arisen from a cause not covered by this limited warranty. This limited warranty does not apply to: (A) Normal routine service items; (B) Repair or replacement necessitated by misuse, abuse, accident, or repairs made by persons other than MVE or persons not authorized by MVE; (C) Defects caused by the effects of normal wear and tear; (D) Acts of God, or other causes not within the control of MVE.

If purchaser believes that a product does not comply with the limited warranty stated above, Purchaser should contact MVE, describing the problem and providing proof of the date of purchase. If directed by MVE Purchaser shall return Products freight prepaid, properly packaged in a MVE approved shipping container and properly identified by a Return Material Authorization Number issued by MVE. Products returned without a Return Material Authorization Number will be refused and returned at Purchaser's expense.

The remedies available for any breach of this limited warranty are limited to repair or replacement of the defective Product or refund of the purchase price at the sole discretion of MVE. MVE warrants that replacement or repaired Products shall be free from defects in material and workmanship for the duration of the unexpired portion of the original warranty or ninety (90) days from the date of re-shipment to Purchaser, whichever is longer. MVE SHALL HAVE NO FURTHER LIABILITY FOR DAMAGES, LOSSES, COST OR FEES OF ANY KIND OR NATURE, WHETHER FORESEEABLE OR NOT, INCLUDING BUT NOT LIMITED TO ATTORNEY'S FEES AND CONSEQUENTIAL, GENERAL, SPECIAL, EXEMPLARY OR PUNITIVE DAMAGES, ARISING OUT OF OR RELATED TO THE USE OF MVE EQUIPMENT, EVEN IF MVE HAS BEEN ADVISED OF THE POSSIBILITY OF SUCH DAMAGES, LOSSES, COST OR FEES.

**EXCEPT FOR THIS LIMITED WARRANTY, MVE HAS MADE NO WARRANTIES OR REPRESENTATIONS, EXPRESSED, OR IMPLIED, INCLUDING BUT NOT LIMITED TO IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. NO REPRESENTATION OR STATEMENT OF MVE MAY CHANGE OR ALTER THIS LIMITED WARRANTY.**

Any claims for breach of this limited warranty shall be governed by Minnesota law and must be brought in a state or federal court in Minnesota.

Some states do not allow limitations on implied warranties or on incidental or consequential damages, so the above limitations may not apply. This limited warranty gives the purchaser specific legal rights. Purchaser may also have other rights, which vary from state to state.

**GENERAL DESCRIPTION**

The SC, XC and LAB series container is a vacuum insulated container of aluminum with fiberglass neck construction providing the highest efficiency possible in liquid nitrogen storage. Use the container for inert fluids only. Liquid oxygen is not compatible with fiberglass material and must not be stored. A sharp blow to the outer vessel can damage the neck tube or start a vacuum leak. Use caution and common sense in handling the container. The life expectancy of the CryoSystem Series is five (5) years. SC, XC, and Lab Series have a life expectancy of ten (10) years.

Upon receipt of the container, examine both the container and packaging for any evidence of damage during shipping. Contact the shipping agent if any damage is apparent. Watch after the first fill for any signs of vacuum loss, such as frost or sweating on the outside jacket. (Some frost near the tip just after filling is normal).

**A. SAFETY**

-  To avoid injury by frostbite, use extreme care whenever handling liquid nitrogen, liquid nitrogen storage, transfer vessels or any objects which have come in contact with liquid nitrogen.
  - Leave no areas of skin exposed.
  - Always wear proper safety attire over clothing: face shield, cryogenic gloves, cryogenic apron.
  - Never overfill liquid nitrogen vessels.
-  Always keep liquid nitrogen vessel in an upright position.
-  Do not tightly seal liquid nitrogen container or prevent nitrogen gas from escaping.
-  Use extreme care to prevent spilling and splashing liquid nitrogen during transfer.
- Immediately remove any clothing or safety attire on which liquid nitrogen has been spilled.
- Get immediate medical attention for any frostbite injuries due to liquid nitrogen.

 **WARNING:** The venting of nitrogen vapors will deplete oxygen in the air, possibly leading to asphyxiation or even death. **DO NOT** store or use liquid container in areas that have poor ventilation.

**B. FILLING INSTRUCTIONS**

- To avoid damage to the aluminum cryogenic vessel which may result in premature vacuum loss, it is important to follow this procedure during the addition of liquid nitrogen to a warm vessel and on subsequent additions.
1. Slowly pour liquid nitrogen into new or warm vessels with a funnel or transfer line when possible.
  2. Allow liquid nitrogen to settle in vessel for 2 hours to completely cool inner.
  3. Fill the vessel to the desired level after the 2 hr settling (cooldown) time.
  4. If filling the dewar from a pressurized source, make sure that the source tank is at a low pressure (1.52 bar or below).
  5. If transfer hose is used for extracting liquid nitrogen from a pressurized liquid source always use a phase separator on the end of the hose.
  6. Remember to always wear proper safety attire over clothing; face shield, cryogenic gloves, and apron.
  7. Never overfill your dewar with liquid nitrogen. Overfilling the tank may cause pumpout components to leak leading to immediate or premature vacuum failure to occur.

**C. MEASURING LIQUID NITROGEN QUANTITY**

1. Use wooden or plastic dipstick. Never use a hollow tube to measure liquid nitrogen.
2. Level will be indicated by frostline, which develops when dipstick is removed.

**D. LIQUID WITHDRAWAL**

1. Liquid withdrawal from LAB units must be accomplished by pouring or utilizing a withdrawal device. Withdrawal device pressurizes to approximately 0.35 bar and the pressure forces liquid up the withdrawal tube out the valve.
2. Always wear proper safety attire; shield, gloves and apron.

**WARRANTY ACTIVATION**

**IMPORTANT! PLEASE FILL OUT THIS REGISTRATION CARD TO ACTIVATE YOUR WARRANTY**

YOUR NAME \_\_\_\_\_  
 ADDRESS \_\_\_\_\_  
 CITY \_\_\_\_\_ STATE \_\_\_\_\_ ZIP \_\_\_\_\_  
 NAME OF DISTRIBUTOR \_\_\_\_\_  
 ADDRESS \_\_\_\_\_  
 SERIAL NUMBER \_\_\_\_\_ DATE PURCHASED \_\_\_\_\_

*What is being stored in the container?*

- Semen                       Embryos                       Skin                       Vaccines                       Blood  
 Bone Marrow                       Cell Lines                       Other

*Was the new container purchased for*

- Replacement                       New Lab                       Project                       Additional Storage

*How did you become aware of MVE products?*

- Advertisement                       Sales Distributor                       Colleague                       Trade Show  
 Thomas Register                       Internet                       Other \_\_\_\_\_

*What Journal/Magazines are you currently receiving?*

(Rate 1 as most useful, 3 as less useful)

1 \_\_\_\_\_ 2 \_\_\_\_\_ 3 \_\_\_\_\_

*How important were the following features that lead to the purchase of this product?*

	Very Important							Not Important		
Quality of freezer	10	9	8	7	6	5	4	3	2	1
MVE reputation	10	9	8	7	6	5	4	3	2	1
Price	10	9	8	7	6	5	4	3	2	1
Storage Capacity	10	9	8	7	6	5	4	3	2	1
Warranty	10	9	8	7	6	5	4	3	2	1
Standard Features	10	9	8	7	6	5	4	3	2	1
Customer Service	10	9	8	7	6	5	4	3	2	1
Distributor Reputation	10	9	8	7	6	5	4	3	2	1
Colleague Referral	10	9	8	7	6	5	4	3	2	1
Temperature Under Lid	10	9	8	7	6	5	4	3	2	1
Nitrogen Consumption	10	9	8	7	6	5	4	3	2	1
Serviceability	10	9	8	7	6	5	4	3	2	1
Other	10	9	8	7	6	5	4	3	2	1

*How satisfied are you with your local MVE representative?*

	Very Satisfied							Not Satisfied		
	10	9	8	7	6	5	4	3	2	1

Comments \_\_\_\_\_

*Overall, how satisfied are you with your MVE container?*

	10	9	8	7	6	5	4	3	2	1

Comments \_\_\_\_\_

Please make any additional comments or suggestions \_\_\_\_\_

**CHART BIO-MEDICAL PRODUCTS**  
**ATTN: CUSTOMER SERVICE MANAGER**  
**112 Windmill Road**  
**Sunbury on Thames**  
**Middlesex**  
**TW167HB UK**

PLACE  
STAMP  
HERE

---

---

---

---

*Thank You!* You have just purchased the finest cryobiological storage container manufactured today. With Chart Bio-Medical Division's experience and reputation, you can be assured of the most advanced technology and unsurpassed quality throughout our product. These containers have been designed, manufactured and tested to achieve maximum efficiency, along with durability to give many years of trouble-free service. If you have any questions regarding this product please contact Chart Bio-Medical.

## WARRANTY

MVE warrants to the original purchaser ("Purchaser") that each new Liquid Nitrogen Dewar (each a "Product"), from the date of shipment, shall be free from defects in materials and workmanship for a period of two (2) years, except as provided below, and dewar vacuum integrity shall be free from defects in materials and workmanship as listed below, except as provided below:

Spectrum and CryoSystem Series, ET11, ET24, A-200, Workhorse 2003: 3 years / XC, SC, and Lab Series, ET 23-JLM: 5 years

MVE warrants all repaired equipment and factory blemished equipment shall be free of defects in materials and workmanship for a period of (90) days from the date of shipment and warrants all dewar vacuum integrity shall be free from defects in materials and workmanship for a period of (1) year from date of shipment. All accessory items are warranted for (90) days from date of shipment to be free of defects in material and workmanship. Purchaser agrees that before this limited warranty shall become effective, Purchaser shall fully inspect each Product within ten (10) days of delivery and before such Product is put to use. Purchaser also agrees to operate the Product in accordance with MVE's operating instructions and that failure to do so shall void this limited warranty. Purchaser further agrees that any claim for breach of warranty must be made in writing within 60 days of discovery of a purported defect. MVE will not be responsible for any alleged breach of warranty, which, as a result of MVE's inspection, MVE determines to have arisen from a cause not covered by this limited warranty. This limited warranty does not apply to: (A) Normal routine service items; (B) Repair or replacement necessitated by misuse, abuse, accident, or repairs made by persons other than MVE or persons not authorized by MVE; (C) Defects caused by the effects of normal wear and tear; (D) Acts of God, or other causes not within the control of MVE.

If purchaser believes that a product does not comply with the limited warranty stated above, Purchaser should contact MVE, describing the problem and providing proof of the date of purchase. If directed by MVE Purchaser shall return Products freight prepaid, properly packaged in a MVE approved shipping container and properly identified by a Return Material Authorization Number issued by MVE. Products returned without a Return Material Authorization Number will be refused and returned at Purchaser's expense.

The remedies available for any breach of this limited warranty are limited to repair or replacement of the defective Product or refund of the purchase price at the sole discretion of MVE. MVE warrants that replacement or repaired Products shall be free from defects in material and workmanship for the duration of the unexpired portion of the original warranty or ninety (90) days from the date of re-shipment to Purchaser, whichever is longer. MVE SHALL HAVE NO FURTHER LIABILITY FOR DAMAGES, LOSSES, COST OR FEES OF ANY KIND OR NATURE, WHETHER FORESEEABLE OR NOT, INCLUDING BUT NOT LIMITED TO ATTORNEY'S FEES AND CONSEQUENTIAL, GENERAL, SPECIAL, EXEMPLARY OR PUNITIVE DAMAGES, ARISING OUT OF OR RELATED TO THE USE OF MVE EQUIPMENT, EVEN IF MVE HAS BEEN ADVISED OF THE POSSIBILITY OF SUCH DAMAGES, LOSSES, COST OR FEES.

**EXCEPT FOR THIS LIMITED WARRANTY, MVE HAS MADE NO WARRANTIES OR REPRESENTATIONS, EXPRESSED, OR IMPLIED, INCLUDING BUT NOT LIMITED TO IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. NO REPRESENTATION OR STATEMENT OF MVE MAY CHANGE OR ALTER THIS LIMITED WARRANTY.**

Any claims for breach of this limited warranty shall be governed by Minnesota law and must be brought in a state or federal court in Minnesota.

Some states do not allow limitations on implied warranties or on incidental or consequential damages, so the above limitations may not apply. This limited warranty gives the purchaser specific legal rights. Purchaser may also have other rights, which vary from state to state.

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Medical Product Services, Borngasse 20, 35619 Braunsfels, Germany

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